Vakuumtechnik im System
Technology for Vacuum Systems

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Instructions for repair: 20999112 - 07 / version 01.07.2018

Valid for Diaphragm pumps
Vacuum gauges (except of VAP 5)

It is the users' responsibility to check the validity of the pump's specific drawing and spare parts list based on the serial number of the pump.

If the pump is older or newer please contact VACUUBRAND and order the pump's specific drawing and spare parts list stating the pump type, serial number and year of manufacture.

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As part of the implementation process of our new ERP system in 2018 our catalog numbers are extended to 8 digits in length by adding the digits “2” or “20” to the beginning of the old catalog numbers.

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☞ Warning! Possible danger. Severe injuries as well as damage to equipment and environment can occur.

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• Caution! Possible danger. Slight injuries as well as damage to equipment and environment can occur.

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Note. Disregarding of notes may cause damage to the product.

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Netzstecker ziehen.
Isolate equipment from mains before removing the cover.

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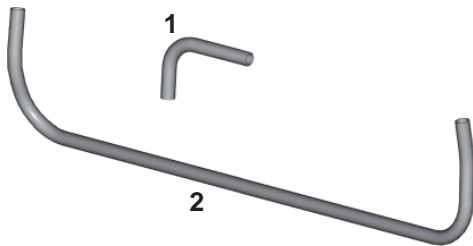
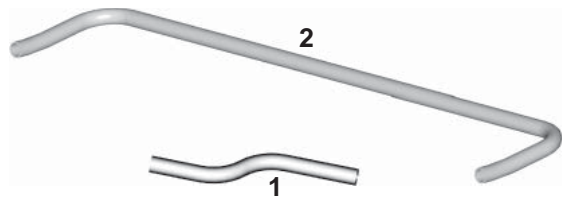
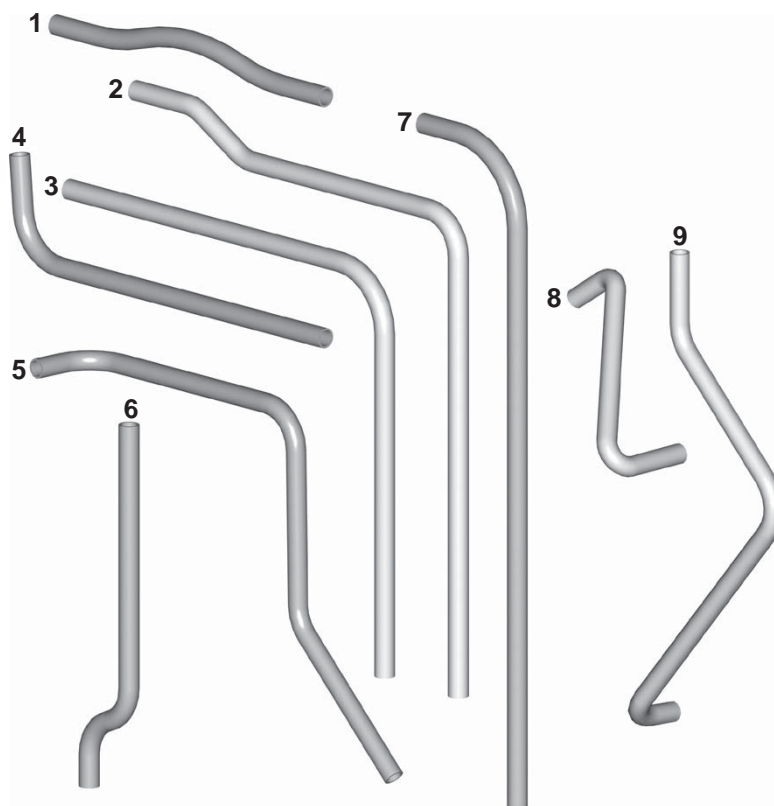
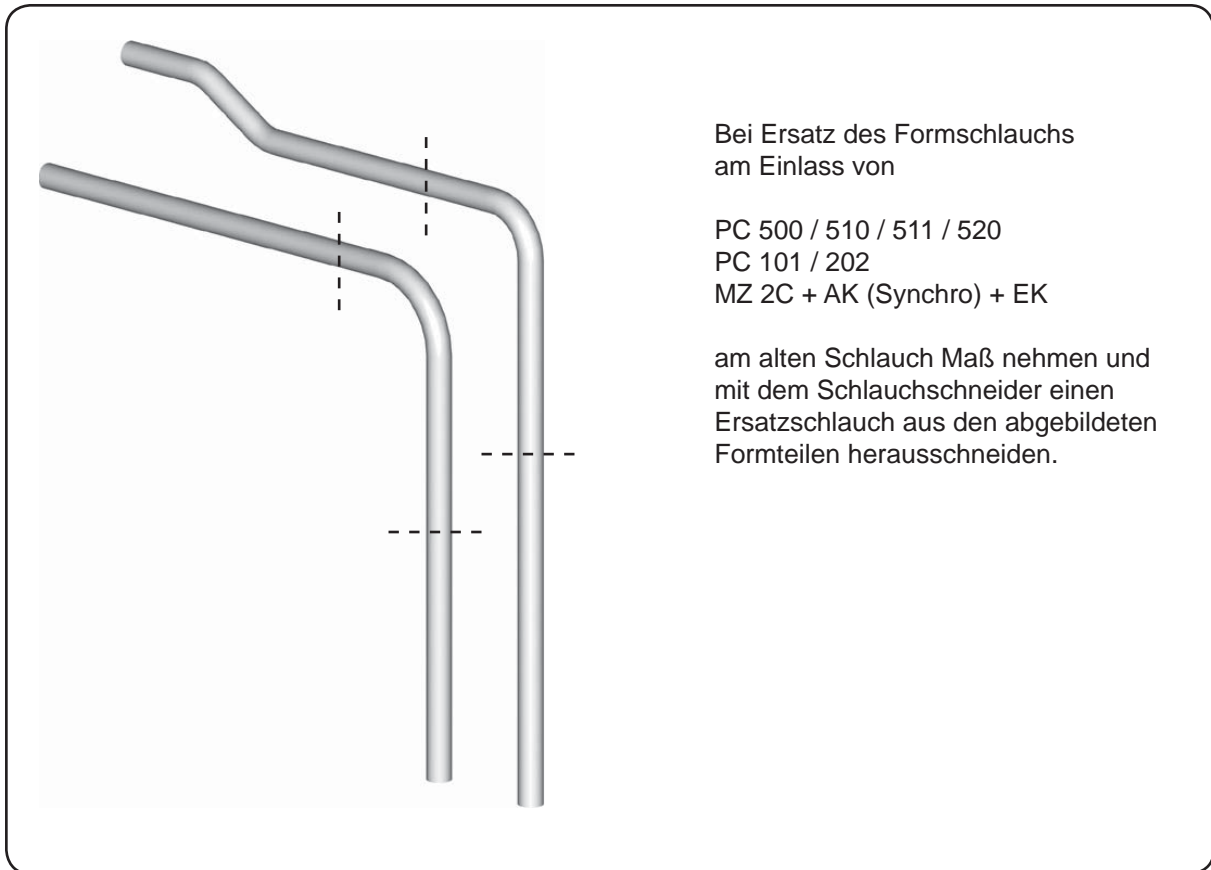
Safety information!

NOTICE

Ensure that repair is done only by an appropriately trained and supervised technician. Obey local and national safety requirements (safety at work law and regulations for environmental protection). Ensure that the technician is familiar with the safety procedures which relate to the products processed by the pumping system. Ensure that the pump is decontaminated before repair and that you take adequate precautions to protect people from the effects of hazardous substances if contamination has occurred.

Read the **instructions for use** of the equipment carefully, especially the sections "Safety information" and "Use and operation". You may order the applicable "Instructions for use" from VACUUBRAND or download at www.vacuubrand.com.

Before starting repair, check, that the required OEM spare parts are available and of the correct type. Check that the required tools are at hand.

► Use exploded view drawings, spare parts lists and circuit diagrams to check the operating sequence notionally beforehand on feasibility, safety requirements and consequences on safety and operability of the equipment.

WARNING

⚠ Before starting repair, isolate the pump and other components from the vacuum system and the electrical supply. Unplug the power cord. Allow the pump to cool down, so that it is at safe temperature for skin contact.

⚠ Before starting maintenance, **wait two minutes** after isolating the equipment from mains to allow the capacitors to discharge.

⚠ Ensure that the pump cannot be operated accidentally. **Never operate the pump if covers or other parts of the pump are disassembled. Never operate a defective or damaged pump.**

CAUTION

If **hazardous or corrosive substances** have been pumped:

- The pump might be contaminated with the process chemicals which have been pumped during operation. Ensure that the pump is decontaminated before maintenance.
- Take adequate precautions to protect people from the effects of dangerous substances. Wear appropriate safety-clothing. Do not inhale and avoid skin contact.
- Obey regulations when disposing of **solvents and chemicals**. Keep in mind that solvents and chemicals might be contaminated with pumped substances.
- Use only **OEM spare parts and accessories**. Otherwise safety and performance of the equipment as well as the electromagnetic compatibility of the equipment might be reduced. Possibly the CE mark or the cTÜVus mark become void if not using OEM spare parts.
- It is mandatory to replace any damaged parts.
- Check operability and safety after repair.
 - If any problems arise, return the pump to the factory for inspection or repair if necessary.

In order to comply with law (occupational, health and safety regulations, safety at work law and regulations for environmental protection) vacuum pumps, components and measuring instruments returned to the manufacturer can be repaired only when certain procedures (see section "**Notes on return to the factory**") are followed.

Scrapping and waste disposal:

Dispose of the pump and any components removed from it safely in accordance with all local and national safety and environmental requirements.

Particular care must be taken with components and waste oil which have been contaminated with dangerous substances from the process.

⚠ You may authorize us to dispose of the equipment **at your expense**.

General notes

Inspection and maintenance of the pump heads and **replacement of diaphragms and valves**: See instructions for use of the diaphragm pump.

All parts of the pump head can be replaced individually, the connecting rod unit is an assembly and must be replaced completely as such.

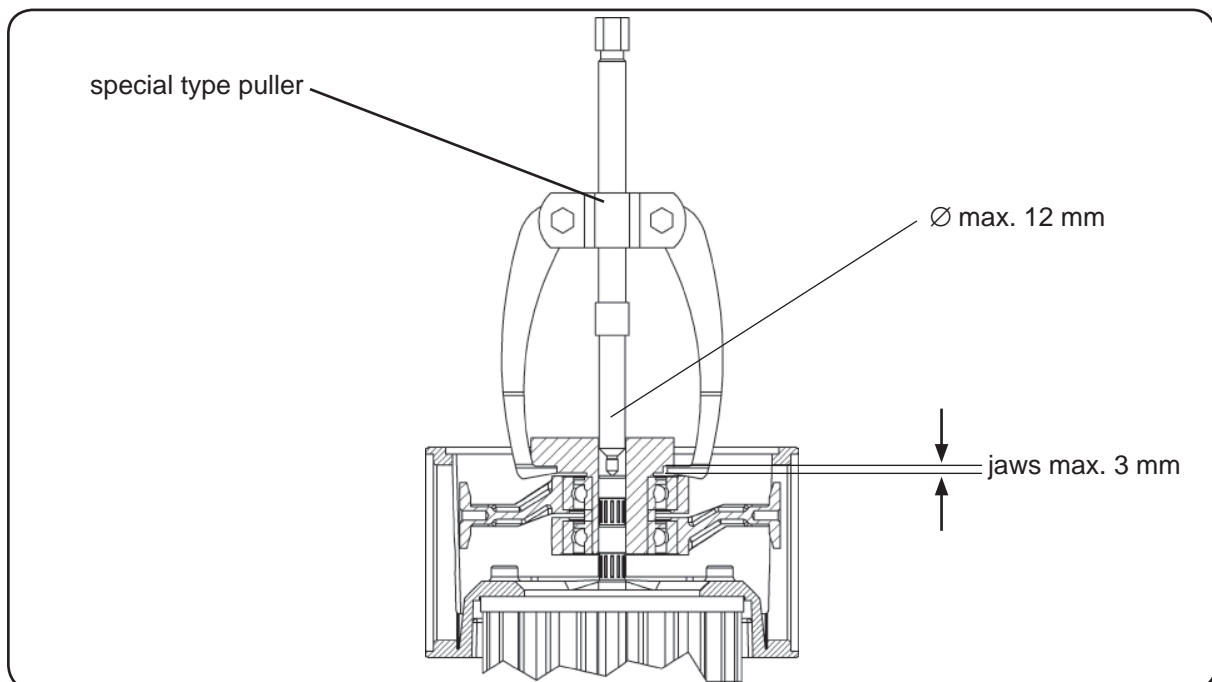
Tools required (details for all types of pumps):

- Hex key size 5
- Open-ended wrench w/f 14, 16, 17
- Phillips screw driver size 2
- Face wrench size 40/4
- **Special type puller** (only for pumps with one, two or four cylinders) with lengthening piece, with jaws max. 3 mm (**cat.-no.: 2613381**)

Replacing the connecting rods

Replacement of the connecting rods of the **ME 16(C)**, **MZ 8(C)**, **MD 8(C/VARIO)**, **MD 12(C)** or **MV 10(C/VARIO)** pumps should be carried out in the factory only.

- ➔ Disassemble the pump heads as described in the instructions for use.
- ➔ Use the special type puller and remove eccentric bushing with connecting rod.
- 📖 Note position of connecting rod relative to housing bore.

- ➔ Assemble the new connecting rods in the reverse order of the disassembling procedure. Note the following:
 - ➔ Mount new tolerance rings on the motor shaft.
 - ➔ Use a suitable press to install the connecting rods (e. g. bench drill machine or lever press machine).
- 📖 Do not exert any force on the motor bearing. Hence the motor has to be supported via the motor shaft while the rods are being installed. Therefore hold up the motor shaft, e.g. with a mandrel, at the center of the motor shaft (recess at end of motor shaft) at the side of the motor fan (pumps with one or two cylinders) or at the opposite end of the motor shaft (pumps with four cylinders; removing of the opposite housing plate of the pump is required).
- 📖 Make sure connecting rods are installed in correct position relative to bore.

- ☞ When press-fitting the connecting rods, make sure that each rod is exactly centred with respect to the corresponding housing bore. Position the rods onto the motor shaft and press in but not completely, press only so far until the diaphragms can be assembled through the housing bores. The rods may still be off-center. Assemble diaphragms with diaphragm support disc and diaphragm clamping disc. Then check the diaphragms' bearing surfaces on the housing. Press in the rods until the diaphragms are centered exactly with respect to housing bores.
- ☞ Pumps with four cylinders: Install the second pair of connecting rods at the opposite side of the pump shifted by 180°.

Attention: If the diaphragm is not correctly centred with respect to the housing bore, the lifetime of the diaphragms and bearings will be reduced and may lead to a leak caused by pulling the diaphragm out of its clamping at the outer diameter.

Replacing the motor

- ➔ Remove the connecting rod as described in section "Replacing the connecting rod".
- ☞ If necessary remove the fan housing of the separately driven fan, pay attention to the cables between fan and terminal box.
- ➔ Remove socket head screws and remove the diaphragm pump housing from the motor.
- ☞ In case, disassemble the terminal box from the defective motor and assemble it on the new motor.
- ➔ Check mains voltage and type of current (see rating plate).
- ➔ Mount the pump housing onto the motor and fix with socket head screws. In case, assemble fan housing of a separately driven fan. Press-fit the connecting rods.

Adjusting the number of washers after replacing the rod or the motor

- ☞ Check performance of the individual pump heads.
- ☞ Allow the pump to warm up (approx. 20 minutes).
- ➔ Measure the pressure at the inlet port of the individual head: Use a suitable, correctly calibrated vacuum gauge (e. g. DVR 2, cat. no.: 682902) and measure the pressure at the inlet port. A vacuum of less than 85 mbar should be indicated.
- ☞ If the reading is higher, increase the number of washers between connecting rod and diaphragm support disc.
- ☞ If the pump at operating temperature becomes noisy, reduce the number of washers.

Notes regarding the use of the sets of hoses

Attention: The sets of hoses are designated for different types of pumps. Therefore the set of hoses consists of hoses, which may not be necessary in the individual case.

Redundant parts are not credited by VACUUBRAND.

The sets of hoses consist of formed hoses and cut goods.

When replacing straight hoses, measure the old hose and cut off the appropriate length using the enclosed hose cutter.

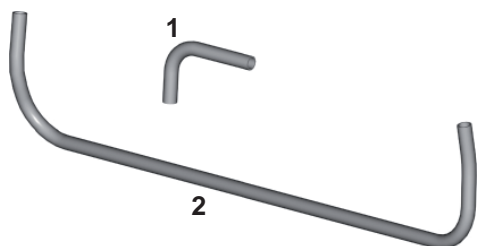
Position union nuts on the hoses. If necessary widen the ends of the hoses prior to assembly by warming (hoses made of PE max. 150°C / hoses made of PTFE max. 280°C).

Attention: The blade of the hose cutter is very sharp. Risk of injury!


⚠ CAUTION

**set of hoses 696850
(only formed hoses)**


- 1: MV 10
MD 12
- 2: ME 16 (2x)
MZ 8
MD 8
MD 12 (2x)
MV 10

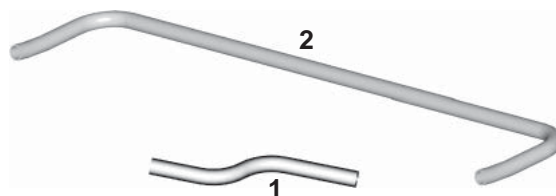**set of hoses 696851
(only formed hoses)**


- 1: ME 16C
MZ 8C
MD 8C
- 2: ME 16C (2x)
MZ 8C
MD 8C
MD 12C (2x)
MV 10C

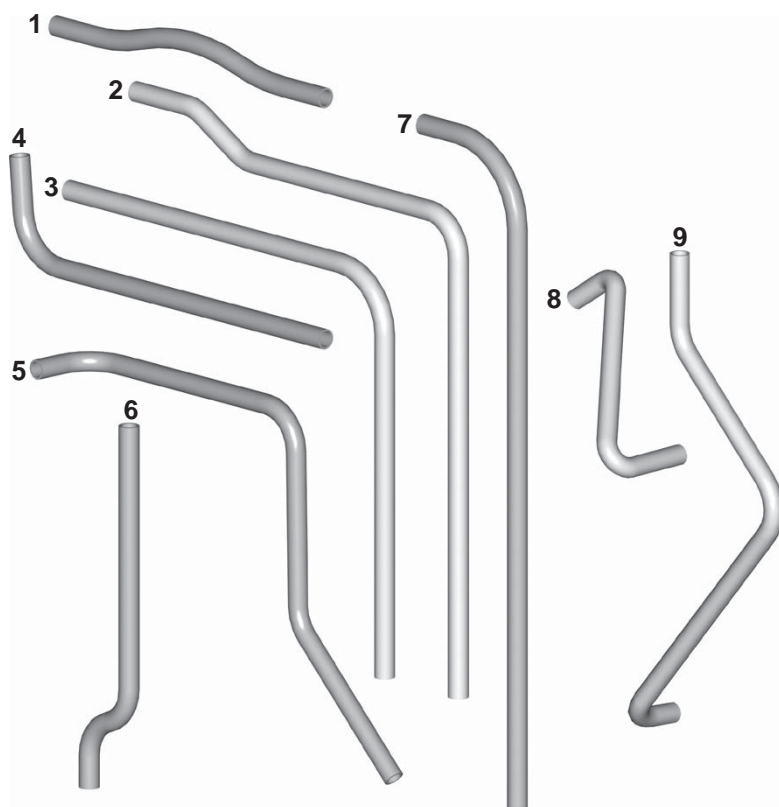**set of hoses 696852
(only formed hoses)**

- 1: PC 2004 VARIO
- 2: PC 2004 VARIO
PC 600 / 610 / 611 / 620
MD 4C + AK (Synchro) + EK
- 3: PC 2002 VARIO
- 4: PC 2003 VARIO
- 5: PC 2003 VARIO
PC 2004 VARIO
PC 600 / 610 / 611 / 620
MD 4C + AK (Synchro) + EK
- 6: PC 600 / 610 / 611 / 620
MD 4C + AK (Synchro) + EK
- 7: PC 2008 VARIO
PC 2010 VARIO
PC 8 / MD 8C
PC 8 / MV 10C
- 8: PC 500 / 510 / 511 / 520
MZ 2C + AK (Synchro) + EK
- 9: PC 2002 VARIO

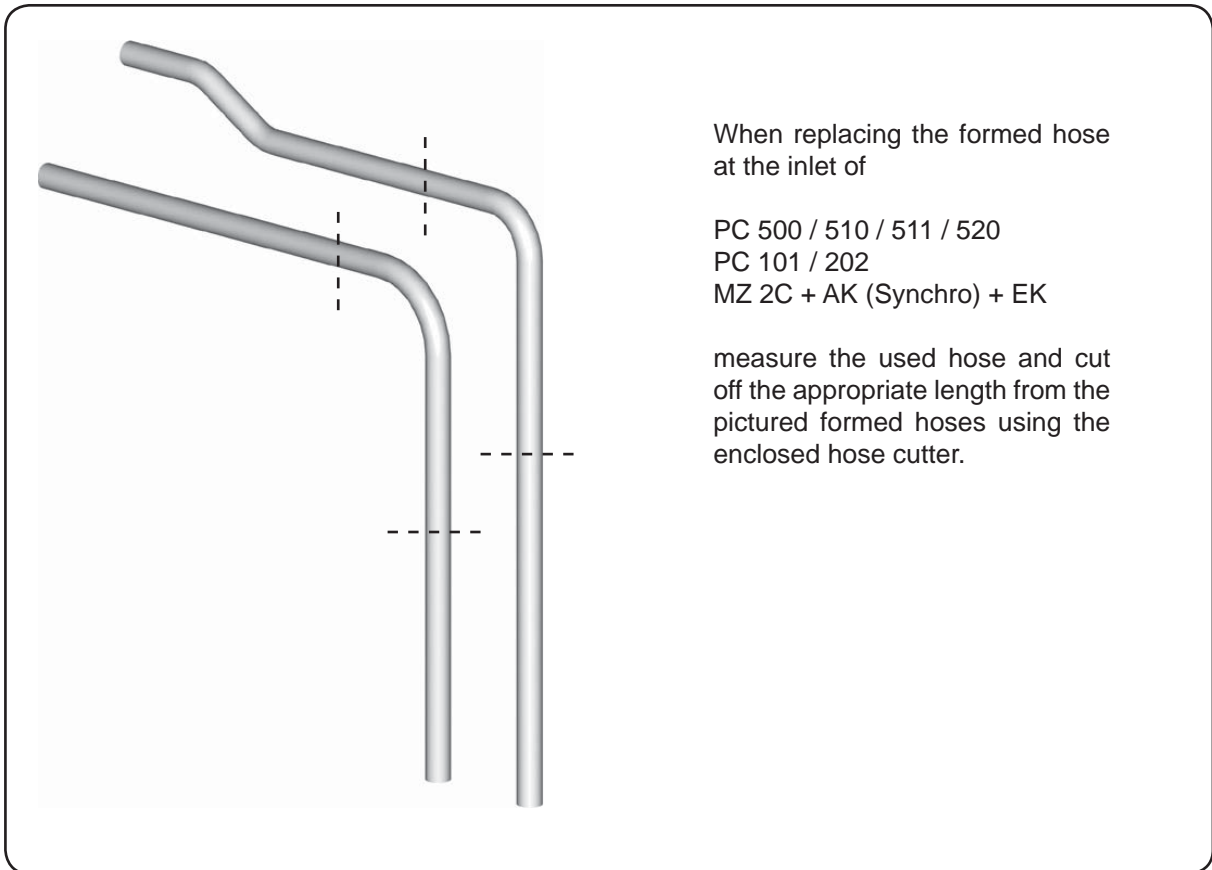When replacing the formed hose
at the inlet of

PC 500 / 510 / 511 / 520

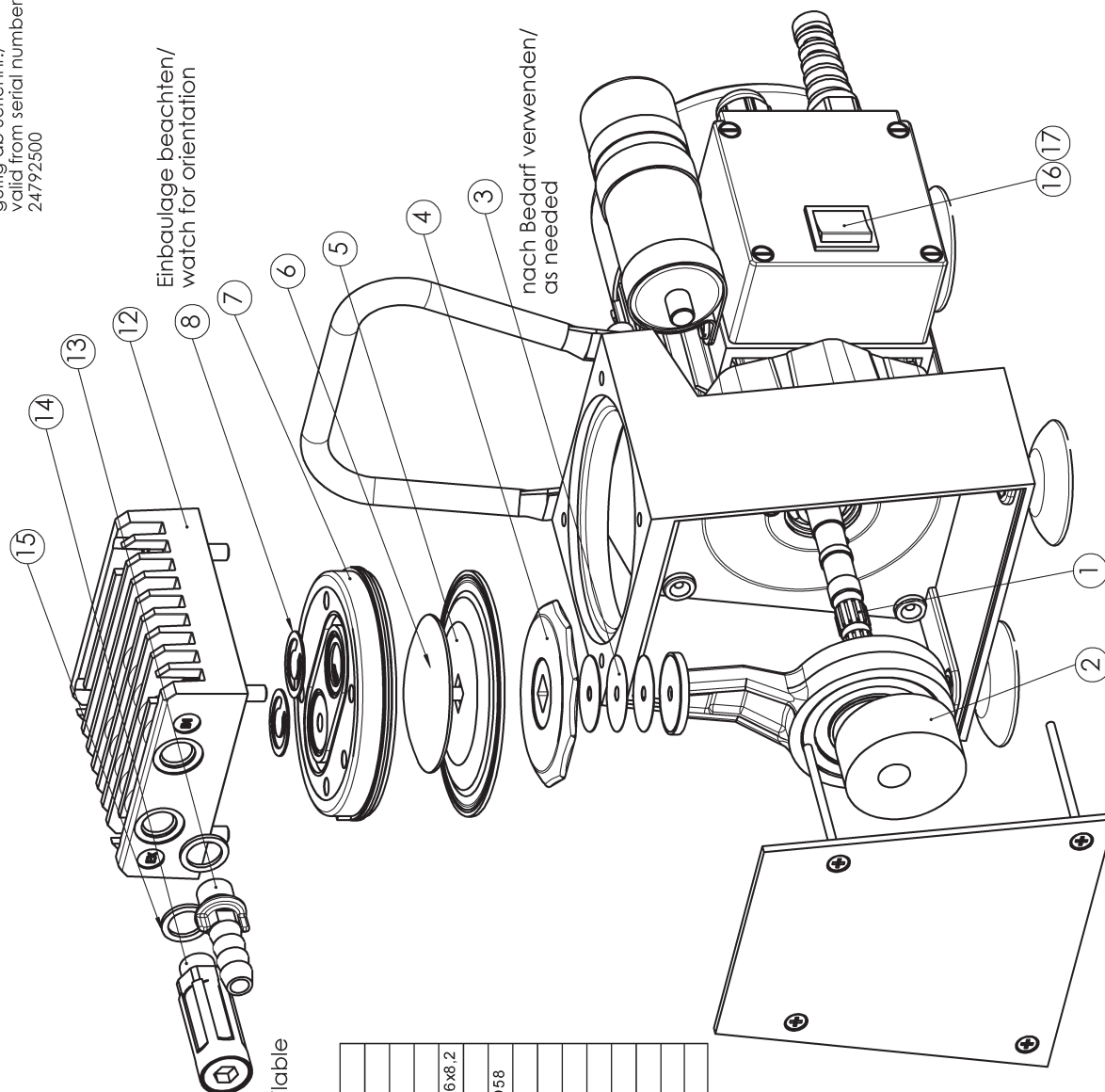
PC 101 / 202

MZ 2C + AK (Synchro) + EK

measure the used hose and cut
off the appropriate length from the
pictured formed hoses using the
enclosed hose cutter.

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DVR 2

Open the equipment

- ➔ Unscrew the four countersunk head screws at the rear side of the housing using a Phillips screw driver.
- ➔ Disassemble the housing parts carefully.

When replacing the batteries be aware to match the positive and negative polarities (as shown by the engraving inside the compartment)!

Replacing the batteries

- ➔ Remove the used batteries / battery and replace by batteries / battery of the recommended type (lithium Mignon-cells (AA), Energizer®* FR 6 1.5V or from serial number 22164301 onwards battery 9V, lithium, 1.2 Ah, Ultralife U9VL). Dispose of old batteries according to regulations.

* reg. trade mark Ralston

Replacing the contact springs

(not applicable from serial number 22164301 onwards)

- ➔ Remove the batteries. Replace the defective contact springs in the battery compartment by new ones.

Replacing the pcb with LCD

(only from serial number 22164301 onwards)

- ➔ Remove the plug of the connecting cable from the pcb. **Attention:** Do not pull at the connecting cable.
- ➔ Replace the defective pcb with LCD by a new one.
- ➔ Attach cable to pcb.

Replacing the LCD

- ➔ Remove the housing front part with glass pane.
- ➔ Unplug the defective LCD from the pcb and plug in the new LCD, make sure that the mark at the LCD is on the left side. When assembling the LCD, only exert pressure to the side with the pins of the LCD.

Replacing the housing front part with glass pane

- ➔ Remove the defective housing front part with glass pane and replace it by a new one.

Replacing the rear side of the housing with connecting flange and pressure transducer

- ➔ Remove the plug of the connecting cable from the pcb. **Attention:** Do not pull at the connecting cable.
- ➔ Attach the connecting cable of the rear side of the housing to the pcb.

Reassembling the equipment

- ☞ When reassembling the vacuum gauge, make sure that the switch pad and the pcb are positioned correctly.
- ➔ Carefully screw both parts of the housing together.

After repair readjust the device (see instructions for use)!

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CVC 2^{II} / CVC 2000 / DVR 5

Before starting repair, **isolate the equipment** from the vacuum system and **the electrical supply**.

Tools required: Phillips screwdriver size 2

How to open the equipment

- ➔ Unscrew the three countersunk screws at the rear side of the housing, the four screws at the bottom of the equipment and both binding screws at the housing front part using a Phillips screw driver.
- ➔ Remove the housing rear part of the equipment (if necessary, remove hose nozzle with O-ring), remove the housing and open the equipment carefully.
- ⚠ Note: The parts on the printed circuit board are sensitive to electrostatic discharge.
- ➔ Do not touch components.
- ➔ If necessary, connect printed circuit board and technician to ground.
- ➔ Remove ribbon contact from the power supply's printed circuit board (pcb) and from the reading's printed circuit board.

Reassembling

- ➔ After having finished the repair, reassemble the ribbon cable and screw the housing together.

After repair, readjust the device (see instructions for use)!

CVC 2^{II}

Replacing the pressure transducer

- ➔ Remove the hose from the pressure transducer.
- ➔ Remove the defective pressure transducer and install the new one.
- ➔ Attach hose to the pressure transducer.

Replacing the pressure transducer hose

- ➔ Remove the hose from the pressure transducer and from the valve flange.
- ➔ Attach new pressure transducer hose.

Replacing the valve flange with valve

- ➔ Remove the pressure transducer hose from the valve flange.
- ➔ Remove the plug of the valve flange from the power supply pcb.
- ➔ Unscrew the two Phillips screws and remove the valve flange.
- ➔ Insert and tighten the new valve flange.
- ➔ Attach the plug of the valve flange to the power supply pcb.
- ➔ Attach the pressure transducer hose to the valve flange.

Replacing the mains filter with switch

- ➔ Remove the three plug-in connections of the mains filter from the power supply pcb.
- ➔ Unscrew the two Phillips screws and remove the mains filter.
- ➔ Assemble the new mains filter.
- ➔ Plug the three plug-in connections of the mains filter onto the power supply pcb.

Replacing the ribbon cable

- ➔ Remove the ribbon cable from the power supply pcb and from the reading pcb.
- ➔ Assemble the new ribbon cable.

Replacing the reading pcb

- ➔ Remove the hose from the pressure transducer.
- ➔ Remove the plugs of both foil cables from the reading pcb.
- ➔ Unscrew the four Phillips screws and remove the defective reading pcb from the housing front part.
- ➔ Screw the new reading pcb with LCD and the housing front part together.
- ➔ Plug in both foil cables.
- ➔ Attach the hose of the pressure transducer.

Replacing the power supply pcb

- ➔ Remove the plugs of the valve flange and of the mains filter from the power supply pcb.
- ➔ Unscrew the Phillips screws and remove the defective power supply pcb from the lower part of the housing.
- ➔ Fasten the new power supply pcb with screws.
- ➔ Attach the plugs of the valve flange and of the mains filter onto the power supply pcb.

Replacing the EPROM

- ➔ Remove the defective EPROM and replace it by a new one.

Replacing the plugged-in LCD

- ➔ Remove the hose from the pressure transducer.
- ➔ Remove the plugs of both foil cables from the reading pcb.
- ➔ Unscrew the four Phillips screws and remove the defective reading pcb from the front side of the housing.
- ➔ Unplug the defective LCD from the pcb and assemble the new LCD, make sure that the mark at the LCD is at the side of the VACUUBRAND logo on the reading pcb. When assembling the LCD, only exert pressure to the side with the pins of the LCD.
- ➔ Screw the reading pcb with LCD and the housing front part together.
- ➔ Attach the foil cables.
- ➔ Attach the hose to the pressure transducer.

Replacing the componentry housing front part with glass pane and reading pcb with plugged-in LCD (for controllers without plugged-in LCD)

- ➔ Remove the hose from the pressure transducer.
- ➔ Attach the hose to the pressure transducer of the pressure transducer of the new reading pcb (assembled to the housing front with glass pane, foil keys and attached LCD).

Replacing the componentry housing front part with glass pane (for controllers with plugged-in LCD)

- ➔ Remove the plugs of both foil cables from the reading pcb.
- ➔ Unscrew the four Phillips screws and remove the reading pcb together with the plugged-in LCD.
- ➔ Assemble the reading pcb with LCD to the new housing front part with glass pane.
- ➔ Attach the foil cables.

CVC 2000 / DVR 5

Replacing the mains filter with switch

- ➔ Remove the three plug-in connections of the mains filter from the power supply pcb.
- ➔ Unscrew the two Phillips screws and remove the mains filter.
- ➔ Assemble the new mains filter.
- ➔ Plug the three plug-in connections of the mains filter onto the power supply pcb.

Replacing the ribbon cable

- ➔ Remove the ribbon cable from the power supply pcb and from the reading pcb.
- ➔ Assemble the new ribbon cable.

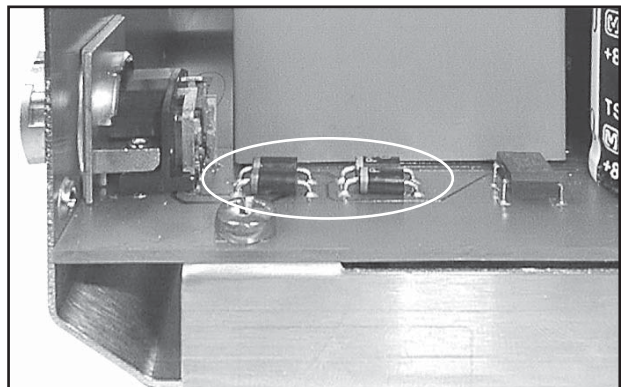
Replacing the read out pcb

- ➔ Unscrew the four Phillips screws and remove the defective reading pcb from the housing front part.
- ➔ Screw the new reading pcb with LCD and the housing front part together.

Replacing the power supply pcb

- ➔ Remove the three plug-in connections of the mains filter from the power supply pcb.
- ➔ Unscrew the Phillips screws (in case, remove nut from the sensor plug; only new versions) and remove the defective power supply pcb from the lower part of the housing.
- ➔ Fasten the new power supply pcb with screws.
- ➔ Plug the three plug-in connections of the mains filter onto the power supply pcb (fasten nut at the sensor plug if available).

Attention, DVR 5 only: Remove four diodes (see fig.) before assembling the new power supply pcb.

Replacing the EPROM

- ➔ Remove the defective EPROM (note correct position) and replace it by a new one.

Replacing the plug-in LCD

- ➔ Unscrew the four Phillips screws and remove the defective reading pcb from the front side of the housing.
- ➔ Unplug the defective LCD from the pcb and assemble the new LCD, make sure that the mark at the LCD is at the side of the VACUUBRAND logo on the reading pcb. When assembling the LCD, only exert pressure to the side with the pins of the LCD.
- ➔ Screw together the reading pcb with LCD and the housing front part.

Replacing the componentry housing front part with glass pane and read out pcb with plug-in LCD (for controllers without plug-in LCD)

- ➔ Remove defective component group by a new one.

Replacing the componentry housing front part with glass pane (for controllers with plug-in LCD)


- ➔ Unscrew the four Phillips screws and remove the reading pcb together with the plug-in LCD.
- ➔ Assemble the reading pcb with LCD to the new housing front part with glass pane.

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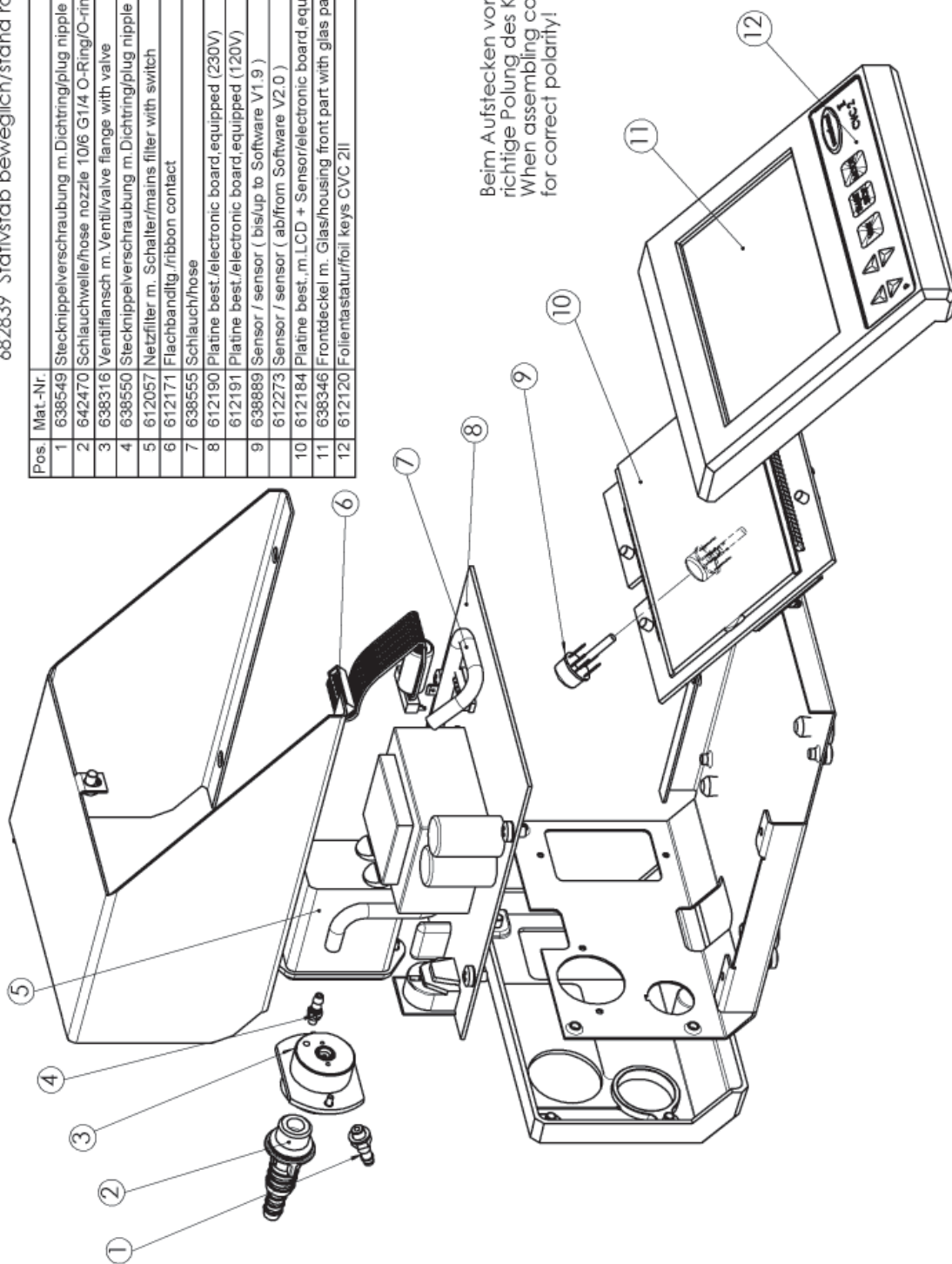Repair - Maintenance - Return - Calibration

IMPORTANT

Every employer (user) is held responsible for the health and safety of his employees. This also applies to service personnel performing repair, maintenance, return or calibration.

The **health and safety clearance form** informs the contractor about any possible contamination of the device and forms the basis for the risk assessment.

In case of devices which have been in contact with biological substances of risk level 2 contact the VACUUBRAND service absolutely before dispatching the device. These devices have to be completely disassembled and decontaminated by the user prior to shipment. Do not return devices which have been in contact biological substances of risk level 3 or 4. These devices cannot be checked, maintained or repaired. Also decontaminated devices must not returned to VACUUBRAND due to a residual risk.

The same conditions apply to on-site work.

No repair, maintenance, return or calibration is possible unless the correctly completed health and safety clearance form is returned. Devices sent are rejected if applicable. Send a completed copy of the **health and safety clearance form** to us in advance. The declaration must arrive before the equipment. Enclose a second completed copy with the product.

Remove all components from the device that are no original VACUUBRAND components. VACUUBRAND will not be responsible for lost or damaged components that are no original components.

Drain the device completely of fluids and residues. Decontaminate the device.

Close all openings airtight especially if using substances hazardous to health.

To expedite repair and to reduce costs, please enclose a detailed description of the problem and the product's operating conditions with every product returned.

If you do not wish a repair on the basis of our **quotation**, the device may be returned to you disassembled and at your expense.

In many cases, the components must be cleaned in the factory prior to repair.

For cleaning we use an environmentally friendly water based process. Unfortunately the combined attack of elevated temperature, cleaning agent, ultrasonic treatment and mechanical stress (from pressurised water) may result in damage to the paint. Please mark in the health and safety clearance form if you wish a repaint at your expense just in case such a damage should occur. We will also replace parts for cosmetic reasons at your request and at your expense.

Before returning the device

Pack the device properly, if necessary, please order original packaging materials at your costs.

Mark the package completely

Enclose the completed health and safety clearance form.

Notify the carrier of any possible contamination if required.

Scrapping and waste disposal

Dispose of the equipment and any components removed from it safely in accordance with all local and national safety and environmental requirements. Particular care must be taken with components and waste oil which have been contaminated with dangerous substances from your processes. Do not incinerate fluoroelastomer seals and O-rings. You may authorize us to dispose of the equipment **at your expense**. Otherwise we return the device at your expense.

Health and safety clearance form

1. Device (Model):
2. Serial no.:
3. Reason for return / malfunction:
4. Has the device been used in a copper process step (e.g., semiconductor production): yes no
5. Substances (gases, liquids, solids, biological material, e. g. bacteria, viruses) in contact with the device / which have been pumped:

6. Risk level of the used biological material: none 1 2* 3** 4**
 * Contact the VACUUBRAND service absolutely before dispatching the device.
 ** Devices which have been in contact with biological substances of risk level 3 or 4 cannot be checked, maintained or repaired. Also decontaminated devices must not returned to VACUUBRAND due to a residual risk.
7. Radioactive contamination: yes no
8. Prior to return to the factory the device has been decontaminated: yes no
 Description of the decontamination method and the test / verification procedure:

9. All parts of the device are free of hazardous, harmful substances: yes no
10. Protective measures required for service staff:

11. If the paint is damaged, we wish a repaint or a replacement of parts for reason of appearance (repaint and replacement at customer's expense): yes no

12. Legally binding declaration

We assure for the returned device that all substances, which have been in contact with the device are listed in section 5 and that the information is complete and that we have not withheld any information. We declare that all measures - where applicable - have been taken listed in section "Repair - Maintenance - Return - Calibration". By our signature below, we acknowledge that we accept liability for any damage caused by providing incomplete or incorrect information and that we shall indemnify VACUUBRAND from any claims as regards damages from third parties. We are aware that as expressed in § 823 BGB (Public Law Code of Germany) we are directly liable for injuries or damages suffered by third parties, particularly VACUUBRAND employees occupied with handling/repairing the product. Shipping of the device must take place according to regulations.

Name: Signature:

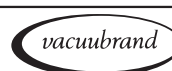
Job title: Company's seal:

Date:

Release for repair grant by VACUUBRAND (date / signature):.....
 Protective measures: Protective gloves, safety goggles Hood External cleaning

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Disclaimer: Our technical literature is only intended to inform our customer. The validity of general empirical values and results obtained under test conditions for specific applications depend on a number of factors beyond our control. It is therefore strictly the users' responsibility to very carefully check the validity of application to their specific requirements. No claims arising from the information provided in this literature will, consequently, be entertained.

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© 2018 VACUUBRAND GMBH + CO KG Printed in Germany

Alfred-Zippe-Str. 4 · 97877 Wertheim / Germany
T +49 9342 808-0 · F +49 9342 808-5555
info@vacuubrand.com · www.vacuubrand.com

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